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ADCP Fc γ RIIb Jurkat Effector Cell Line

ADCP Fc γ RIIb Jurkat Effector 细胞系

For research use only!

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Reagent	Specification	Manufacturer/Catalogue No.
Puromycin	25 mg	Genomeditech/GM-040401-1
Blasticidin	10 mg	Genomeditech/GM-040404-1
Pen/Strep(P.S)	100 mL	Thermo/15140-122
Fetal Bovine Serum	500 mL	Cegrogen biotech/A0500-3010
RPMI 1640	500 mL	Viva Cell BIOSCIENCES/C3010-0500
96 Well Clear V-Bottom Tissue Culture	96-well	Corning/3894
96 well round well culture plate	96-well	NEST/701001
96 well White Flat Bottom Polystyrene Not Treated Microplate	96-well	Corning/3912
Anti-H_MS4A1(CD20) hIgG1 Antibody(Ocrelizumab)	/	Genomeditech/GM-27200AB
APC anti-human CD32B/C	100 tests	Biologend/398304
Raji Cell Line	/	Genomeditech/GM-C19100
GMOne-Step Luciferase Reporter Gene Assay Kit	1000T	Genomeditech/GM-040503C

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Equipment	Manufacturer/Catalogue No.
细胞计数仪	ThermoFisher Scientific/Countess 3
酶标仪	Moleculardevices/SpectraMax L

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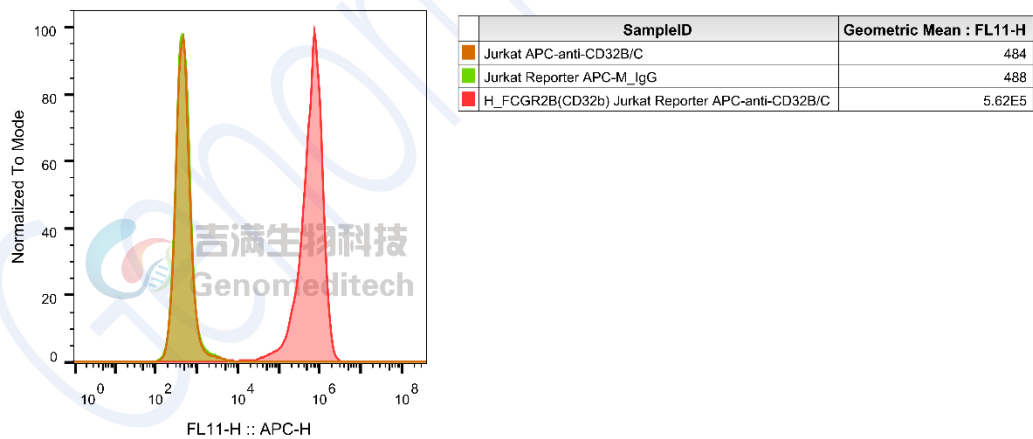
Fig 4. 使用 APC anti-human CD32B/C (Biolegend/398304) 流式验证结果

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FcγR	
Cynomolgus_FcRn MDCK Cell Line	H_FCGR1A(CD64) CHO-K1 Cell Line
H_FCGR1A(CD64) HEK-293 Cell Line	H_FCGR2A(CD32A) CHO-K1 Cell Line
H_FCGR2B(CD32B) CHO-K1 Cell Line	H_FCGR3A(CD16a) 158F CHO-K1 Cell Line
H_FCGR3A(CD16a) 158V CHO-K1 Cell Line	H_FCGR3B(CD16b) CHO-K1 Cell Line
H_FcRn CHO-K1 Cell Line	H_FcRn MDCK Cell Line
Mouse_FcRn MDCK Cell Line	
Anti-FcRn hIgG4 Reference Antibody(Rozabio)	Anti-H_FcRn IgG4 Antibody(Rozanolixizumab)
Anti-Mouse CD1632 mIgG2b Antibody(2.4G2)	
ADCCP	
ADCC FcγRIIIa(158F) Jurkat Effector Cell Line	ADCC FcγRIIIa(158V) DDX35TM Jurkat Effector Cell Line
ADCC FcγRIIIa(158V) Jurkat Effector Cell Line	ADCC M_FcγRIV Jurkat Effector Cell Line
ADCP FcγRIIa DDX35TM Jurkat Effector Cell Line	ADCP FcγRIIa Jurkat Effector Cell Line
ADCP FcγRIIa R131 Jurkat Effector Cell Line	

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