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ADCP Fc γ RI Jurkat Effector Cell Line

ADCP Fc γ RI Jurkat Effector 细胞系

For research use only!

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Reagent	Specification	Manufacturer/Catalogue No.
Puromycin	25 mg	Genomeditech/GM-040401-1
Blasticidin	10 mg	Genomeditech/GM-040404
Pen/Strep(P.S)	100 mL	Thermo/15140-122
Fetal Bovine Serum	500 mL	Cegrogen biotech/A0500-3010
RPMI 1640	500 mL	Viva Cell BIOSCIENCES/C3010-0500
96 Well Clear V-Bottom Tissue Culture	96-well	Corning/3894
96 well round well culture plate	96-well	NEST/701001
96 well White Flat Bottom Polystyrene Not Treated Microplate	96-well	Corning/3912
Anti-H ₄ MS4A1(CD20) hIgG1 Antibody(Ocrelizumab)	/	Genomeditech/GM-27200AB
Raji Cell Line	/	Genomeditech/GM-C19100
PE anti-human CD64 Antibody	25 tests	Biologend/305007
GMOne-Step Luciferase Reporter Gene Assay Kit	1000T	Genomeditech/GM-040503C

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Equipment	Manufacturer/Catalogue No.
细胞计数仪	ThermoFisher Scientific/Countess 3
酶标仪	Moleculardevices/SpectraMax L

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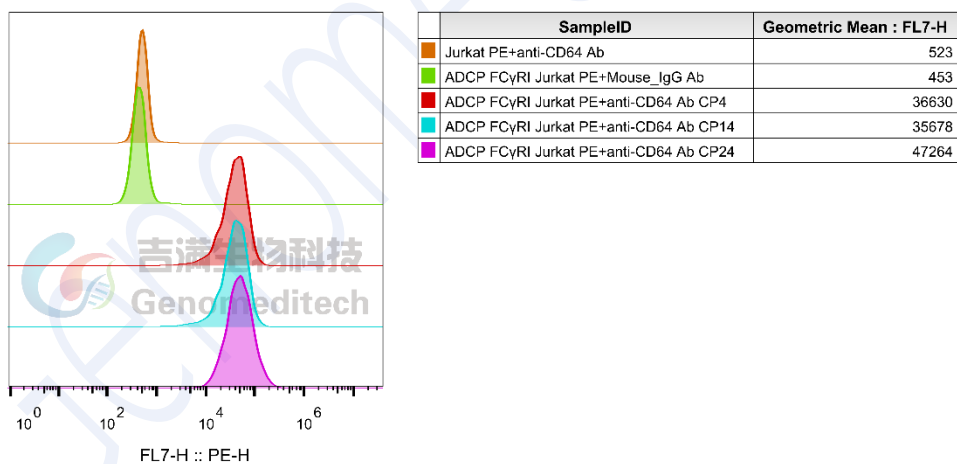
Fig 4. 使用 PE anti-human CD64 Antibody (Biolegend/305007)抗体流式稳定性验证结果

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FcγR	
Cynomolgus_FcRn MDCK Cell Line	H_FCGR1A(CD64) CHO-K1 Cell Line
H_FCGR1A(CD64) HEK-293 Cell Line	H_FCGR2A(CD32A) CHO-K1 Cell Line
H_FCGR2B(CD32B) CHO-K1 Cell Line	H_FCGR3A(CD16a) 158F CHO-K1 Cell Line
H_FCGR3A(CD16a) 158V CHO-K1 Cell Line	H_FCGR3B(CD16b) CHO-K1 Cell Line
H_FcRn CHO-K1 Cell Line	H_FcRn MDCK Cell Line
Mouse_FcRn MDCK Cell Line	
Anti-FcRn hIgG4 Reference Antibody(Rozabio)	Anti-H_FcRn IgG4 Antibody(Rozanolixizumab)
Anti-Mouse CD1632 mIgG2b Antibody(2.4G2)	
ADCCP	
ADCC FcγRIIIa(158F) Jurkat Effector Cell Line	ADCC FcγRIIIa(158V) DDX35TM Jurkat Effector Cell Line
ADCC FcγRIIIa(158V) Jurkat Effector Cell Line	ADCC M_FcγRIV Jurkat Effector Cell Line
ADCP FcγRIIa DDX35TM Jurkat Effector Cell Line	ADCP FcγRIIa Jurkat Effector Cell Line
ADCP FcγRIIa R131 Jurkat Effector Cell Line	ADCP FcγRIIb Jurkat Effector Cell Line

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